<u>armfield</u>

PCT SERIES: PROCESS CONTROL TECHNOLOGY



Essentials of Process Control – PCT60–64

The Modern Process Control system from Armfield, allows students to investigate the principles of industrial process control, using independent Temperature, Pressure, Flow, Level and Servo/ Pendulum systems.

An advanced digital controller/logger with USB, Wifi, Bluetooth and LAN communications is built into each set of equipment.

Three sets of Windows PC software are supplied free of charge: Basic control, On/Off control and PID control. Software includes all settings and functions and on-screen graphing of Set point, Process value and Kp, Ki, and Kd terms in the Control Equation which is displayed dynamically. Data can be viewed on the internal graph or saved to a CSV file for later analysis.

mfield

PCT60 - Level Control / PCT61 - Flow Control / PCT62 - Temperature Control PCT63 - Pressure Control / PCT64 - Servo Pendulum

> PROCESS CONTROL PID





- Understanding the drive/sensors
- On/Off control systems
- System time constant
- P, PI and PID control
- Zeigler Nichols algorithm

Integral wind up

Derivative filter

PROCESS

mfield

Manual tuning

lssue: 2

- Interfacing with MATLAB/LabVIEW
- Servo pendulum angle control

armfield

UK office - email: sales@armfield.co.uk tel: +44 (0) 1425 478781 (for ROW) USA office - email: info@armfieldinc.com tel: +1 (609) 208-2800 (USA only)

URL: http://www.armfield.co.uk/pct50

We reserve the right to amend these specifications without prior notice. E&OE \odot 2024 Armfield Ltd. All Rights Reserved

armfield.co.uk

Applications

IP

ChE ME

PROCESS

rmfie

PCT SERIES: PROCESS CONTROL TECHNOLOGY



Level control is a fundamental aspect of process control in various industries, including chemical, petrochemical, food and beverage, water treatment, and more. Understanding level control is essential for students pursuing careers in engineering, process automation, or any field related to these industries.

The Armfield PCT60 level control system introduces students to fundamental control concepts, such as feedback control, proportional-integral-derivative (PID) controllers, and closed-loop control systems.

The Level system consists of a reservoir water tank, a variable speed pump, a pressure-based level sensor, and a clear process vessel with a scale. A proportional valve provides the process vessel drain.

An overflow pipe in the process vessel prevents it from being overfilled and the system allows students to adjust the pump speed and valve opening.

Features/Benefits

- ▶ USB, WiFi, Bluetooth & LAN communications supplied as standard
- Supplied software includes Basic control, On/Off control and PID control
- Software dynamically displays Set point, Process value and Kp, Ki, and Kd.
- MATLAB and Labview compatible
- Supplied with full set of manuals and teaching material

Experimental content


- Understanding how to control driving devices
- PID controller
- Zeigler Nichols algorithm
- Understanding the sensors
 On/Off control systems
- Integral wind up
 Derivative filter
- System time constant
- P controllerPI controller
- Manual tuning
- Interfacing with MATLAB/LabVIEW

Ordering specifications

PCT60 Level Control Process

- A level control process trainer, comprising:
- ▶ 4mm x 1.5mm wall clear PVC tube (300mm)
- ▶ 24V PSU 60w 2.5a
- Level process control system assembly
- ▶ Barley pot for process control
- ► USB lead
- Manual Control software allowing low level access to the Drive and Load in each system allowing calibration of sensors and drive systems
- On/Off Control software allowing control of each system with a simple On/Off algorithm, view software based oscillations and to explore the effects of hysteresis
- PID Control software allowing uses to enter values for Kp, Ki, Kd and see how the system reacts to in-putted values

UK office - email: sales@armfield.co.uk tel: +44 (0) 1425 478781 (for ROW) USA office - email: info@armfieldinc.com tel: +1 (609) 208-2800 (USA only)

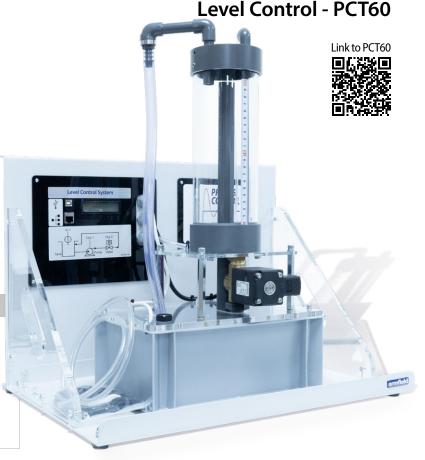



. . . .

Mains electrical supply: 110-230V, AC 50-60 Hz.

PC and Display meeting the following minimum specification:

Scale

- Processor: 1Ghz or faster
- RAM: 1Gb or more
- HDD Space: 1Gb
- **OS:** 32 or 64bit Windows 7, 8, 10 or 11
- Display: Recommended minimum (1920 by 1080) full HD

Overall dimensions

| Length | 64cm | |
|---|---------------------|--|
| Width | 45cm | |
| Height | 33cm | |
| Packed and crated shipping specifications | | |
| Volume | 0.095m ² | |
| Gross weight | 12.1kg | |

Ordering code

PCT60-UK: Level Process Control System PCT60-EU: Level Process Control System PCT60-USA: Level Process Control System

We reserve the right to amend these specifications without prior notice. E&OE © 2024 Armfield Ltd. All Rights Reserved

PCT SERIES: PROCESS CONTROL TECHNOLOGY





Flow process control is a critical aspect of various industries, including chemical, petrochemical, manufacturing, and food production.

Students who understand flow control principles are better prepared for careers in these fields.

Proper flow control is essential for optimizing production processes. Students learn how to regulate the flow of materials or fluids in a way that minimizes waste, maximizes product quality, and improves overall process efficiency.

The PCT61 Flow control system introduces students to fundamental control concepts, such as feedback control, proportional-integralderivative (PID) controllers, and closed-loop control systems.

The Flow system consists of a water tank, variable speed pump, a turbine type flow sensor, an electrically operated proportional valve and a variable area flow meter (rotameter).

This allows students to adjust the flow rate via the pump speed and the valve opening to develop a PID based control system.

Features/Benefits

- ▶ USB, WiFi, Bluetooth and LAN communications supplied as standard
- Supplied software includes Basic control, On/Off control and PID control.
- Software dynamically displays Set point, Process value and Kp, Ki, and Kd
- MATLAB and Labview compatible
- Supplied with full set of manuals and teaching material

Experimental content

- Understanding how to control driving devices
- ▶ PID controller
- Zeigler Nichols algorithm
- Understanding the sensors
- On/Off control systems
- System time constant
- P controller
- PI controller

- Integral wind up
- Derivative filter
- Manual tuning
- Interfacing with MATLAB LabVIEW

Ordering specifications

PCT61 Flow Control Process

- A flow control process trainer, comprising:
- 4mm x 1.5mm Wall clear PVC tube (300mm)
- ▶ 24V PSU 60w 2.5a
- Flow process control assembly
- ► Barley pot for process control
- USB lead
- Manual Control software allowing low level access to the Drive and Load in each system allowing calibration of sensors and drive systems
- On / Off Control software allowing control of each system with a simple on off algorithm, view software based oscillations and to explore the effects of hysteresis
- ▶ PID Control software allowing uses to enter values for Kp, Ki, Kd and see how the system reacts to in-putted values

UK office - email: sales@armfield.co.uk tel: +44 (0) 1425 478781 (for ROW) USA office - email: info@armfieldinc.com tel: +1 (609) 208-2800 (USA only)



Related products

PCT60: Level Process Control System

- PCT62: Temperature Process Control System
- PCT63: Pressure Process Control System
- PCT64: Servo Pendulum Process Control System

| Requirements | | |
|--------------|-----|----------|
| РС | USB | . |

Mains electrical supply:

110-230V, AC 50-60 Hz.

PC and Display meeting the following minimum specification:

Scale

- Processor: 1Ghz or faster
- RAM: 1Gb or more
- HDD Space: 1Gb
- OS: 32 or 64bit Windows 7, 8, 10 or 11
- Display: Recommended minimum (1920 by 1080) full HD

Overall dimensions

| Length | 64cm | |
|---|---------------------|--|
| Width | 45cm | |
| Height | 33cm | |
| Packed and crated shipping specifications | | |
| Volume | 0.095m ² | |
| Gross weight | 12.1kg | |

Ordering code

PCT61-UK: Flow Control System PCT61-EU: Flow Control System PCT61-USA: Flow Control System

We reserve the right to amend these specifications without prior notice. E&OE © 2024 Armfield Ltd. All Rights Reserved

PCT SERIES: PROCESS CONTROL TECHNOLOGY



Temperature Control - PCT62

Temperature control is a critical aspect of many industries, including chemical, pharmaceutical, food and beverage, manufacturing, and HVAC (heating, ventilation, and air conditioning). Understanding temperature control is crucial for students pursuing careers in these industries.

Proper temperature control can help industries reduce energy consumption and operating costs. Students learn how to optimize temperature control systems to minimize energy waste and environmental impact.

The Armfield PCT62 temperature control system introduces students to fundamental control concepts, such as feedback control, proportional-integral-derivative (PID) controllers, and closed-loop control systems.

The temperature process control system includes a heated plate within a duct and a thermocouple. A fan at one end of the duct blows ambient air over the block, to change the control conditions and provide a disturbance to the system.

The system allows users to adjust the heater power and the air flow rate to develop a PID based control system then adjust these parameters to achieve the required time/temperature change profilefor the system in response to step changes in system requirements.

Features/Benefits

- ▶ USB, WiFi, Bluetooth and LAN communications supplied as standard
- Supplied software includes Basic control, On/Off control and PID control
- Software dynamically displays Set point, Process value and Kp, Ki, and Kd

PID controller

Integral wind up

Derivative filter

Manual tuning

I abVIFW

Zeigler Nichols algorithm

Interfacing with MATLAB/

- MATLAB and Labview compatible
- Supplied with full set of manuals and teaching material

Experimental content

- Understanding how to control driving devices
- Understanding the sensors
- On/Off control systems
- ► System time constant
- ► P controller
- ► PI controller


Ordering specifications

PCT62 Temperature Control Process

A temperature control process trainer, comprising:

- 24V PSU 60w 2.5a
- Pressure process control assembly
- USB lead
- Manual Control software allowing low level access to the Drive and Load in each system allowing calibration of sensors and drive systems
- On/Off Control software allowing control of each system with a simple On/Off algorithm, view software based oscillations and to explore the effects of hysteresis
- PID Control software allowing uses to enter values for Kp, Ki, Kd and see how the system reacts to in-putted values

UK office - email: sales@armfield.co.uk tel: +44 (0) 1425 478781 (for ROW) USA office - email: info@armfieldinc.com tel: +1 (609) 208-2800 (USA only)

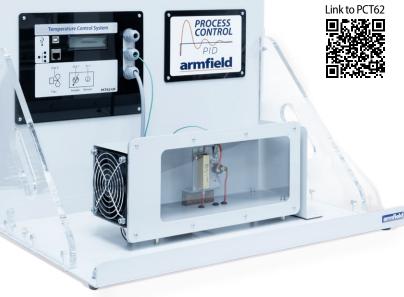

Related products

PCT60: Level Process Control System

- PCT61: Flow Process Control System
- PCT63: Pressure Process Control System
- PCT64: Servo Pendulum Process Control System

| Requirements | Scale |
|--------------------------|-------|
| PC USB | ¢ Î |
| Mains alactrical supply: | |

Mains electrical supply: 110-230V, AC 50-60 Hz.

PC and Display meeting the following minimum specification:

- Processor: 1Ghz or faster
- RAM: 1Gb or more
- HDD Space: 1Gb
- OS: 32 or 64bit Windows 7, 8, 10 or 11
- Display: Recommended minimum (1920 by 1080) full HD

Overall dimensions

| Length | 64cm | |
|---|---------------------|--|
| Width | 45cm | |
| Height | 33cm | |
| Packed and crated shipping specifications | | |
| Volume | 0.095m ² | |
| Gross weight | 12.1kg | |

Ordering code

PCT62-UK: Temperature Control System PCT62-EU: Temperature Control System PCT62-USA: Temperature Control System

We reserve the right to amend these specifications without prior notice. E&OE © 2024 Armfield Ltd. All Rights Reserved

PCT SERIES: PROCESS CONTROL TECHNOLOGY

PCT series

Pressure Process - PCT63

Teaching pressure process control to students equips them with the knowledge and skills necessary for success in industries where pressure regulation is critical.

It provides a strong foundation in engineering, automation, safety, and problem-solving, making students valuable assets in the workforce while contributing to the efficiency, safety, and quality of industrial processes.

Improper pressure levels can lead to equipment failure, leaks, or even explosions and emphasizes the importance of maintaining safe operating conditions.

The Armfield PCT63 system introduces students to fundamental control concepts, as in feedback control, proportional-integral-derivative (PID) controllers, and closed-loop control systems. It consists of a variable-speed reciprocating air pump (compressor), the speed of which can be adjusted by the user, a pressure vessel, and an outflow system.

The outflow system allows the air to escape via either a manually operated needle valve - providing an ongoing outflow, or via a solenoid valve and second needle valve - providing a step change in the outflow.

The vessel pressure is measured by a Bourdon-type mechanical gauge along with a pressure sensor. Providing a visual indication of the pressure in the vessel and a means for the user to check and calibrate the controller input from the pressure sensor.

Features/Benefits

- ▶ USB, WiFi, Bluetooth and LAN communications supplied as standard
- Supplied software includes Basic control, On/Off control and PID control
- Software dynamically displays Set point, Process value and Kp, Ki, and Kd
- MATLAB and Labview compatible
- Supplied with full set of manuals and teaching material

Experimental content

- Understanding how to control driving devices
- Understanding the sensors
- On/Off control systems
- System time constant
- ▶ P controller
- ► PI controller

- ► PID controller
- Zeigler Nichols algorithm
- Integral wind up
- Derivative filter
- Manual tuning
- Interfacing with MATLAB/ LabVIEW

Ordering specifications

PCT63 Pressure Control Process


- A Pressure control process trainer, comprising:
- > 24V PSU 60w 2.5a
- ► Pressure process control assembly
- ► USB lead
- Manual Control software allowing low level access to the Drive and Load in each system allowing calibration of sensors and drive systems
- On/Off Control software allowing control of each system with a simple On/Off algorithm, view software based oscillations and to explore the effects of hysteresis
- ▶ PID Control software allowing uses to enter values for Kp, Ki, Kd and see how the system reacts to in-putted values

UK office - email: sales@armfield.co.uk tel: +44 (0) 1425 478781 (for ROW) USA office - email: info@armfieldinc.com tel: +1 (609) 208-2800 (USA only)



Related products

PCT60: Level Process Control System PCT61: Flow Process Control System PCT62: Pressure Process Control System PCT64: Servo Pendulum Process Control System

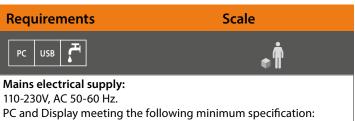

- Processor: 1Ghz or faster
- RAM: 1Gb or more
- HDD Space: 1Gb
- OS: 32 or 64bit Windows 7, 8, 10 or 11
- Display: Recommended minimum (1920 by 1080) full HD

Overall dimensions

| Length | 64cm | |
|---|---------------------|--|
| Width | 45cm | |
| Height | 33cm | |
| Packed and crated shipping specifications | | |
| Volume | 0.095m ² | |
| Gross weight | 12.1kg | |

Ordering code

PCT63-UK: Pressure Process Control System PCT63-EU: Pressure Process Control System PCT63-USA: Pressure Process Control System

We reserve the right to amend these specifications without prior notice. E&OE © 2024 Armfield Ltd. All Rights Reserved

<u>armfield</u>

PCT SERIES: PROCESS CONTROL TECHNOLOGY



Servo Pendulum - PCT64

A servo pendulum is a specific type of control system used for applications like pendulum control and inverted pendulum control.


Learning about servo pendulum control systems allows students to see real-world applications of control theory. These systems are used in areas like robotics, aviation, and mechatronics, and understanding them can help students relate theory to practice.

The Armfield PCT64 servo pendulum control system introduces students to fundamental control concepts, such as feedback control, proportional-integral-derivative (PID) controllers, and closed-loop control systems.

The system consists of a powerful DC motor mounted on a rugged frame. A disc with captive nuts is attached to the DC motor and students are able to screw in 100-gram weights to different parts

of the disc to alter the characteristics of the system.

A single weight at 0 degrees forms an inverting pendulum.

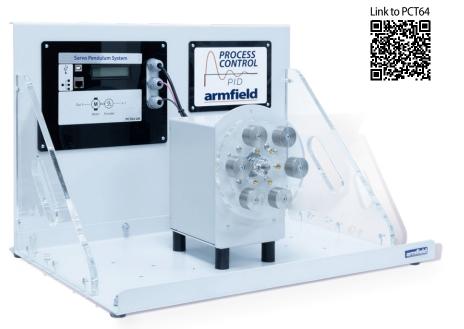

Features/Benefits

- ▶ USB, WiFi, Bluetooth and LAN communications supplied as standard
- Supplied software includes Basic control, On/Off control and PID control
- Software dynamically displays set point, Process value and Kp, Ki, and Kd
- MATLAB and Labview compatible
- Supplied with full set of manuals and teaching material

Experimental content

- Understanding how to control driving devices
- Understanding the sensors
- On/Off control systems
- System time constant
- P controller
- PI controller

- PID controller
- Zeigler Nichols algorithm
- Integral wind up
- Derivative filter
- Manual tuning
- Interfacing with MATLAB/ LabVIEW

Ordering specifications

PCT64 Servo Process Control System Process

- A Servo Pendulum Process Control trainer, comprising:
- 24V PSU 60w 2.5a
- Servo Pendulum control assembly
- ► USB lead
- Manual Control software allowing low level access to the Drive and Load in each system allowing calibration of sensors and drive systems.
- On/Off Control software allowing control of each system with a simple On/Off algorithm, view software based oscillations and to explore the effects of hysteresis
- ▶ PID Control software allowing uses to enter values for Kp, Ki, Kd and see how the system reacts to in-putted values

UK office - email: sales@armfield.co.uk tel: +44 (0) 1425 478781 (for ROW) USA office - email: info@armfieldinc.com tel: +1 (609) 208-2800 (USA only)

Related products

PCT60: Level Process Control System PCT61: Flow Process Control System PCT62: Temperature Process Control System

- PCT63: Pressure Process Control System



110-230V, AC 50-60 Hz.

PC and Display meeting the following minimum specification:

- Processor: 1Ghz or faster
- RAM: 1Gb or more
- HDD Space: 1Gb
- OS: 32 or 64bit Windows 7, 8, 10 or 11
- Display: Recommended minimum (1920 by 1080) full HD

Overall dimensions

| Length | 64cm | |
|---|---------------------|--|
| Width | 45cm | |
| Height | 33cm | |
| Packed and crated shipping specifications | | |
| Volume | 0.095m ² | |
| Gross weight | 12.1kg | |

Ordering code

PCT64-UK: Servo Pendulum Control System PCT64-EU: Servo Pendulum Control System PCT64-USA: Servo Pendulum Control System

We reserve the right to amend these specifications without prior notice. E&OE \otimes 2024 Armfield Ltd. All Rights Reserved


PCT SERIES: PROCESS CONTROL TECHNOLOGY




SOFTWARE AND INTERFACING FOR THE ESSENTIALS OF PROCESS CONTROL UNITS

Each process is supplied complete with software that allows it to be controlled using a Windows PC via a USB connection. There are currently 3 software applications available with the unit.

Manual Control gives low level access to the Drive and Load in each system and allows students to calibrate the sensors and drive systems.

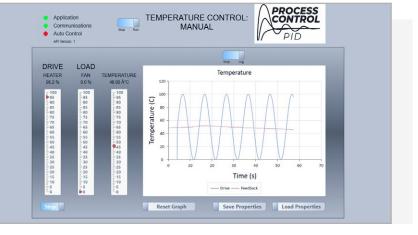

On/Off Control allows students to explore controlling each system with a simple on off algorithm, to see the oscillations this produces and to explore the effects of hysteresis.

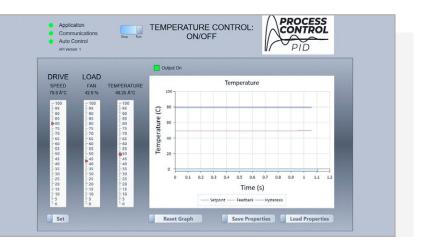

TEMPERATURE CONTROL: PROCESS Communio PID Auto Control PID PID Control allows students to enter values DRIVE LOAD for Kp, Ki, Kd and see how the system reacts Temperature SETPOIN 70.0 Å*C FAN 60.7 % IPERATUR 46.75 ŰC to those values. Control value contributions -100 -95 -90 -85 -80 -75 -50 -55 -50 -45 -30 -25 -20 -15 -10 80 95 90 85 80 75 65 65 50 45 30 25 20 15 10 5 0 95 90 85 80 75 70 65 60 55 50 45 40 35 30 25 20 15 (C) nperature 60 40 -50 20 14 10 12 18 Time (s) Cont 1-0 Set Save Properties Load Properties Reset Graph ERROR 23.25 P 140 I + D 100 -0 Output 100.0 %

The MATLAB software is supplied with the unit. The Control functionality is controlled via a USB COM port or Networked connection to the computer. We have provided a library of functions to allow the various aspects of the Control Hardware to be investigated and controlled. An Application Programming Interface is provided which makes the equipment compatible with MATLAB, Labview or other software.


UK office - email: sales@armfield.co.uk tel: +44 (0) 1425 478781 (for ROW) USA office - email: info@armfieldinc.com tel: +1 (609) 208-2800 (USA only)

We reserve the right to amend these specifications without prior notice. E&OE © 2024 Armfield Ltd. All Rights Reserved

Our Commitment to you

Armfield recognises that it is not enough to just supply quality engineering equipment, but that it must also ensure a complete range of services both pre and post-sale:

- Supplied equipment meets global curriculum requirements
- Expert consultation in laboratory design and layout
- Professional installation and commissioning service
- Comprehensive training for all products in house or on site
- In house trials (industrial and research)
- Two year warranty on all products
- Dedicated aftersales service
- Detailed learning outcomes and experiments supplied with all equipment

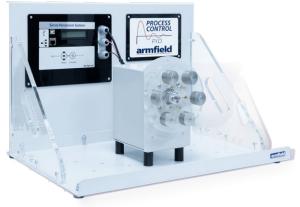

armfield worldwide

REPRESENTATIVES IN MORE THAN 80 COUNTRIES

To locate a contact in your area visit: www.armfield.co.uk/contact-armfield/

U.S. Office:

Armfield Inc. 9 Trenton Lakewood Road Clarksburg NJ 08510 USA

Tel/Fax: +1 (609) 208-2800 E-mail: info@armfieldinc.com Support: armfieldassist.com

Head Office:

Armfield Limited 10 Headlands Business Park, Ringwood, Hampshire BH24 3PB England

Telephone: +44 (0)1425 478781 E-mail: sales@armfield.co.uk Support: armfieldassist.com

Aftercare Installation

Installation Commissioning Training Service and maintenance Support: armfieldassist.com

Knowledge base

> 28 years' expertise in research & development technology
 > 50 years' providing engaging engineering teaching equipment
 Benefit from our experience, just call or email to discuss your
 laboratory needs, latest project or application.

An ISO 9001:2015 Company