


This benchtop unit allows a variety of experiments to be undertaken to investigate test specimens under torsional loading and bending loading within their elastic limits. An extruded base frame carries two solid vertical supports. Profiled grooves within the base frame, fasteners and quick release handles ensure quick and easy adjustment of the span between the vertical supports.

ALLOWS STUDENTS TO INVESTIGATE TEST SPECIMENS UNDER TORSIONAL LOADING AND BENDING LOADING WITHIN THEIR ELASTIC LIMITS

SV807 Dial gauge, 0.01mm resolution

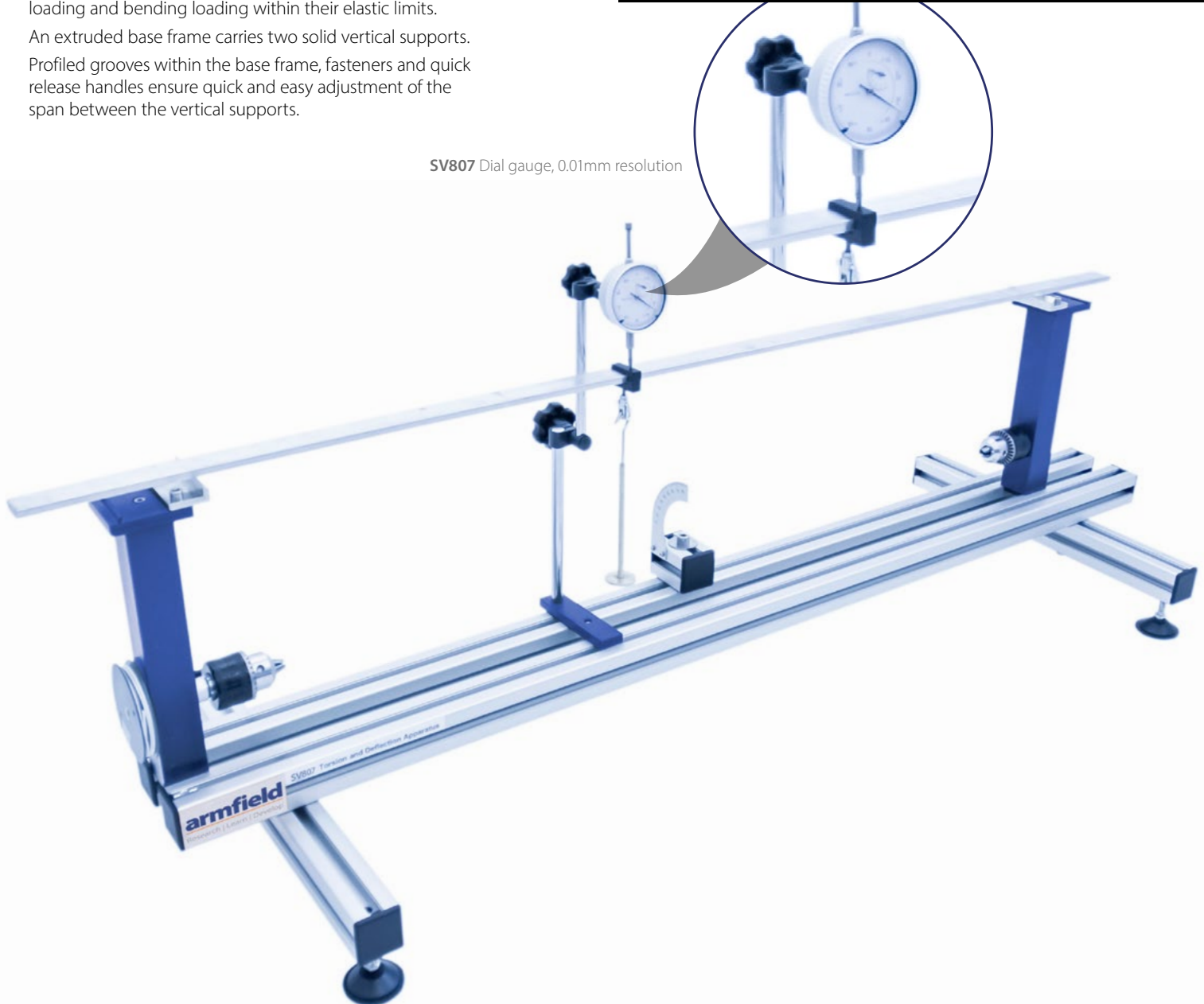

Experimental content

- ▶ Deflection of specimen as a function of loading force, material, Young's Modulus of Elasticity (E), cross section and support span
- ▶ Comparison of bending stiffness of varying specimen sections for the same cross sectional area
- ▶ Comparison of cantilever beams, simply supported beams
- ▶ Determination of Young's Modulus of Elasticity (E) in shear for the different material specimens
- ▶ Introducing Poisson's Ratio
- ▶ Torsion angle and clamping length relationship
- ▶ Torsion angle and torsion moment relationship
- ▶ Torsion angle and specimen cross section
- ▶ Modulus of rigidity introduction

UK office - email: sales@armfield.co.uk tel: +44 (0) 1425 478781 (for ROW)
USA office - email: info@armfield.inc tel: +1 (609) 208-2800 (USA only)

Issue: 1

URL: <http://www.armfield.co.uk/structures>

Applications

ME CE IP

We reserve the right to amend these specifications without prior notice. E&OE © 2022 Armfield Ltd. All Rights Reserved

Description

This benchtop unit allows a variety of experiments to be undertaken to investigate test specimens under torsional loading and bending loading within their elastic limits. An extruded base frame carries two solid vertical supports. Profiled grooves within the base frame, fasteners and quick release handles ensure quick and easy adjustment of the span between the vertical supports.

One vertical support carries a unique specimen chuck with a through hole. Another chuck is held stationary at one end of the base frame.

The chucks have two functions; to secure the torsion specimens during testing and to transmit the torsional loading to the torsion specimen. The torsional load is transmitted using a pulley and cord arrangement, with the pulley being mounted to a shaft running in bearings.

The bearings minimise the friction in the system. The chuck installed into the vertical support has a through hole to allow the length of the torsion specimen to be varied.

The twist of the torsion specimens is indicated on an angle indicator and pointer, which can be moved along the test specimen. The remaining vertical support is used for bending tests.

Each vertical support has a top clamp provided.

This allows the end support arrangement for the bending specimens to be varied, i.e. clamped, simply supported on knife edges or on ground dowels. The supports can be moved along the base frame thus changing the beam span.

To measure the bending specimen deflection a dial gauge is fixed onto a movable stand and can be positioned anywhere within the beam span. A set of torsion and bending specimens are supplied. A load hanger and set of calibrated weights create specimen loading.

Technical specification

- ▶ Torsion specimens: $\text{Ø}5.00 \pm 0.5\text{mm}$ x 470mm long, steel, aluminium, brass and nylon
- ▶ Bending specimens: 1200mm long, steel, aluminium and brass
- ▶ Bending specimens:
 - 25.4mm (W) x 3.18mm (H)
 - 25.4mm (W) x 4.76mm (H)
 - 25.4mm (W) x 6.35mm (H)
 - 19.05mm (W) x 3.18mm (H)
- ▶ 2 x Load hangers
- ▶ 1 x Dial gauge: 25mm travel, 0.01mm resolution, with precision anvil
- ▶ 1 x Movable stand for dial gauge
- ▶ 1 x Tape measure
- ▶ 2 x Clamp plates for cantilevers
- ▶ 2 x Pointer and protractor
- ▶ Knife edge and ground dowels for simply supported beam
- ▶ Torsion shaft running in bearings

Overall dimensions

Length	1.000m
Width	0.243m
Height	0.200m
Packed and crated shipping specifications	
Volume	0.39m ³
Gross weight	53kg

Related laws

- ▶ Mechanical Engineering
- ▶ Structural Engineering
- ▶ Stress
- ▶ Bending
- ▶ Torsion
- ▶ Deflection
- ▶ Span
- ▶ Bending Theory
- ▶ Modulus of Rigidity
- ▶ Super-position

Requirements

Scale



- ▶ Sturdy benchtop

Features / benefits

- ▶ Demonstrates torsion and bending
- ▶ Cantilevers can be tested
- ▶ Specimen bending, deflection
- ▶ Specimen torsion
- ▶ Angle of twist along specimen length
- ▶ Customer specific specimens can be used
- ▶ Three bending specimens
- ▶ Four torsional specimens
- ▶ Full set of weights and hangers supplied

Ordering specification

- ▶ 1 x SV807 Torsion and Deflection Testing Apparatus
- ▶ 4 x Torsion specimen
- ▶ 8 x Bending specimen
- ▶ 2 x Clamp plate
- ▶ 2 x Chuck key
- ▶ 2 x Beam clamp
- ▶ 2 x Load hanger
- ▶ 1 x Hex wrench set
- ▶ 1 x Tape measure
- ▶ 1 x Dial gauge
- ▶ Weights: 5 x 0.2N; 1 x 1N; 2 x 2N; 1 x 5N
- ▶ 1 x 2m Spare cord
- ▶ 1 x Pointer assembly
- ▶ Instruction manual
- ▶ Packing list
- ▶ Test sheet

Operational conditions

- ▶ Storage temperature: -10°C to +70°C
- ▶ Operating temperature range: +10°C to +50°C
- ▶ Operating relative humidity range: 0 to 95%, non-condensing

Ordering codes

- ▶ **SV807:** Torsion and Deflection Testing Apparatus

Armfield standard warranty applies with this product

Knowledge base

- > 28 years expertise in research & development technology
- > 50 years providing engaging engineering teaching equipment

Benefit from our experience, just call or email to discuss your laboratory needs, latest project or application.

An ISO 9001:2015 Company



armfield.co.uk

Aftercare

Installation
Commissioning
Training
Service and maintenance
Support: armfieldassist.com