<u>armfield</u>

Friction

The apparatus consists of a wall mounted pulley with a loaded belt. The pulley is made of aluminium and has two machined grooves to suit a flat, vee belt and a rope. Each belt fits into its respective groove during testing.

For a given belt tension and angle of lap, a turning moment can be applied by adding weights to the pulley drive hanger. The idea is to determine the torque which just causes the pulley to turn, and so to find the ratio of the belt tensions. This enables the belt tension equation to be verified and the coefficient of friction to be determined. The angle of lap can be easily varied by placing one end of each belt at different angular positions on the wall bracket. A removable pin and bar are used to enable the various angular positions to be achieved.

Experimental content

- To observe the relationship between the tensions in the two sides of a belt
- To assess the differences between vee, rope and flat belts
 To determine the coefficient of friction between the pulley
- and belt for the belt sections
- ► To investigate the effect of angle of lap

Related laws

- Automotive
- Torque
- Friction
- Machines
- Wear

Slip

Belt drives

Angle of lap

- hines
- ır

Features / benefits

- Can be mounted to the SV100 Bench Mounted Frame
- One pulley caters for three belts. No interchanging of pulleys
- Constant tangential drive of pulley
- Three belts supplied; flat, vee and rope
- Customer specific belts can be used
- Full set of weights and hanger supplied

Technical specification

- ► Single large pulley
- ► Flat groove and 40° 'V' groove
- ► 150mm effective groove diameter
- Pulley timing grooves at 18° intervals
- Belt anchor points at 15° intervals
- Safety interlock to stop pulley rotating

Overall dimensions

Length	0.230m	
Width	0.150m	
Height	0.280m	
Packed and crated shipping specifications		
Volume	0.01m ³	
Gross weight	20kg	

UK office - email: sales@armfield.co.uk tel: +44 (0) 1425 478781 (for ROW) USA office - email: info@armfield.inc tel: +1 (609) 208-2800 (USA only)

Theory of Machines - SV series

Friction of Belts Apparatus – SV906

WALL OR FRAME MOUNTED APPARATUS TO OBSERVE THE RELATIONSHIP BETWEEN THE TENSIONS IN THE TWO SIDES OF A BELT

Scale

Requirements

►

armfield

Rigid vertical support

Ordering specification

- SV906 Friction of Belts Apparatus
- ▶ 1 x Flat belt
- ► 1 x 'V' belt
- 2 x Hook
- ► 1 x Stopwatch
- ▶ 1 x 5m Spare cord
- ▶ 1 x Load hanger
- ▶ 4 × 0.1N; 3 × 0.2N; 1 × 1N; 2 × 2N; 1 × 5N; 2 × 10N; 4 × 20N; 4 × 50N
- Instruction manual

Recommended accessories / equipment

SV100 Bench Mounted Frame

Ordering codes

- **SV906** Friction of Belts Apparatus
- **SV100** Bench Mounted Frame

Issue: 1 Application			ations
URL: http://www.armfield.co.uk/structures	ME	CE	IP
We reserve the right to amend these specifications without prior notice. E&OE © 2022 Arm	field Ltd.	All Rights	Reserved

armfield.co.uk